APPARATUS

CODE	PART NAME	REMARKS
ATG-00-013	atg-8l	Tag Closed Cup Flash Point Tester, Metal Bath Model, External Chiller excluded
ATG-00-014	atg-8w	Tag Closed Cup Flash Point Tester, Water Bath Model

STANDARD ACCESSORIES

CODE	PART NAME	Qty	REMARKS
ATG-02-411	Test Cup for Tag	1	
210-00-061	O-Ring, G-45	1	For Test Cup
320-00-035	Thermofuse	1	Pack of 5, with Insulating Tube.
	Gas Hose φ9 x φ16, 1.5m, with Hose Bands	1	
ABL-01-401	Windscreen for abl-8/atg-8	1	
	Braided Hose with Wire bands, 1.5m	1	For atg-8w only
	Drain Hose with Hose Band	1	For atg-8w only
ATG-04-021	Flow Control Valve	1	For atg-8w only
210-02-191	Insulated Tubing, 2m	2	For atg-8I only
170-01-051	Hose Band, SPN-8	4	For atg-8I only
MPC-02-032	AC Power Cord		3.0m, For 100 to 120V
MPC-02-033			2.5m, For 220 to 240V
	atg-8w/l Manuals	1	Manuals for atg-8w and atg-8l.

OPTIONAL ACCESSORIES

CODE	PART NAME	REMARKS
070-00-068	Printer, BS2-80TS	Prints out test data and instrument settings.
070-00-281	Mini USB Keyboard	Sample ID entering
	CHILLER	For atg-8l only, Temperature range +5 to 55°C

SUGGESTED SPARES for 2 YEARS

CODE	PART NAME	Qty	CODE	PART NAME	Qty
ATG-02-411	Test Cup for Tag	1	ABL-01-003	Electric Igniter, EL-8	2
210-00-061	O-Ring, G-45	1	ATG-04-031	Temperature Sensor for atg-8	1
320-00-035	Thermofuse	1	ABL-01-012	Flash Detector for abl-8/atg-8	1

We reserve the right of changes without prior notice. Printed in Japan 1203(E)

DISTRIBUTED BY

7-10-3 Ayase, Adachi-ku, Tokyo 120-0005 JAPAN

TEL: +81-3-3620-1711 FAX: +81-3-3620-1713 URL: http://www.tanaka-sci.com

E-mail: overseas-group@tanaka-sci.com

TANA Automatic Petroleum Tester

AUTOMATED TAG CLOSED CUP FLASH POINT TESTER



Model atg-8 series is the automated Tag Closed Cup flash point tester from Tanaka. Based on 45 years experience of manufacturing the automated flash point testers, innovative 8th generation series has been released. The theme of 8th generation is "Ergonomic, Versatile and Premium".

- ERGONOMIC COLOR-COORDINATED DISPLAY FOR HIGH VISIBILITY FOR EVERYBODY (UNIVERSAL DESIGN) AND INTUITIVE SWITCH OPERATION.
- ENHANCED VERSATILITY INCLUDING USB PORT FOR OPTIONAL FLASH MEMORY OR KEYBOARD, PASSWORD PROTECTION AND DATA STRAGE.
- ALL-IN-ONE CHASSIS DESIGN: COMPACT, LIGHT-WEIGHT AND PREMIUM EXTERIOR DESIGN.
- CONFORMS TO STANDARD TEST METHOD (ASTM D56, IP304)

SINGLE ACTION SETTING

The cup cover is permanently mounted on a swing-arm assembly, which allows an easy handling of the test cup cover. It makes free from handling the hot cup cover after completing a test.

The swing arm assembly can be lifted up to vertical position for easier test cup cleaning.

EASY OPERATION

Select a test mode and enter the expected flash point; while the instrument executes the test, you are free to do other lab work. The instrument follows the exact procedures prescribed in the test method, and the completion of the test cycle is signaled by beeps. The test result is brightly shown on the color LCD module.

INTERCHANGEABLE IGNITION SOURCE

atg-8w/atg-8l is equipped with both gas and electric igniter as the ignition source. It only takes a few minutes to switch from gas to electric or vice versa.

DRAFT PROTECTION SHIELD COVER

A FM4910 conforming cover insures the flame integrity. This material shall meet the numerical criterion of each index. (1)FPI(Fire Propagation Index)≤6 (2)SDI(Smoke Damage Index)≤0.4

DATA HANDLING

The unit is equipped to store up to 200 successive test results and export to an optional printer or to a LIMS through an RS-232 port or USB memory.

USB PORT

A USB port allows connecting USB keyboard for easy entering sample name and connecting USB flash memory for data copy to a PC or test conditions/parameters copy to the tester.



COMPACT DESIGN & ENERGY EFFICIENT

atg-8l uses Peltier Cells for sample cooling/heating. Just a small chiller with water suffices the cooling requirement. No methanol is required.

A control part and a test part are put into one body which culminated in achieved in the light weighting of the unit. A liquid-cooling metal bath model and conventional water bath model are available by different measurement temperature range.

SPECIFICATIONS

TYPE	All –in one, microprocessor controlled	
TEST STANDARDS	ASTM D56	
MEASURING RANGE	atg-8w: Ambient to +95°C atg-8l: -20°C to +70°C at cooling medium of 10°C 20°C to +95°C at cooling medium of 50°C	
MEASUREMENT MODES	ASTM CRM, ASTM Normal, ASTM SPE (search), ASTM Rapid and User custom	
ВАТН	atg-8w:Water atg-8l:Metal Block	
SAMPLE COOLING/HEATING	atg-8w:by Plate Heater/by by replacing Bath Liquid atg-8l: by Peltier Devices	
DISPLAY	5.7 inch (117 x 88mm) color LCD with LED back light	
TEMPERATURE SENSOR	PT-100 in stainless steel sheath	
FLASH DETECTOR	CRC Thermocouple	
IGNITION SOURCE	Gas ignition with automatic lighting or electric ignition. Interchangeable	
BAROMETRIC CORRECTION	Automatic correction by built-in barometric pressure sensor	
I/O PORT	RS-232C = 1 channel (for PC or Optional Printer) USB=1 channel (for USB keyboard or USB Flash Memory)	
SAFETY SHUTDOWN	Automatically shuts off and the problem is reported by buzzer and display, in case: (a) EFP+10 °C or 95 °C is reached, (b) temperature sensor is found defective, (c) flash detector is found defective, (d) thermofuse is blown, (e) electric Igniter is blown, (f) built-in battery is found drained out, (g) test cover is not set in place, (h)control computer runs away (no display)	
GAS SUPPLY (When Gas Ignition is used)	LP gas or natural gas (Max. pressure:10kPa)	
POWER SUPPLY	100/120VAC or 220/240 VAC, 50/60Hz	
POWER CONSUMPTION	atg-8w: 500VA MAX. atg-8l: 250VA MAX.	
ELECTRICITY CONSUMPTION	atg-8w: 40Wh for 1 test (about 30 min), 15.12gCO ₂ (@ 0.378kgCO ₂ /1kWh) atg-8l: 30Wh for 1 test (about 30 min), 11.34gCO ₂ (@ 0.378kgCO ₂ /1kWh)	
AMBIENT TEMP. RANGE	10°C to 35°C	
DIMENSIONS & WEIGHT	230mm (W) × 470mm (D) × 390mm (H) 17kg	

CHILLER REQUIREMENTS for atg-8I

TYPE	Circulator with heating/cooling capacities. Open bath type.
TEMPERATURE RANGE	5°C to +55°C or wider
PUMP CAPACITY Pressure:0.2bar minimum Flow rate:2L/min minimum	
COOLING CAPACITY	250W@10°C minimum
HEATER CAPACITY 500W minimum	
BATH VOLUME	3L minimum

Julabo F25-ED is suitable for this specification.